


Warming Platform User's Guide

Time

The default time is one (1) hour, shown as 1.0. You can change the time by using the up/down buttons on the right side of the Warming Platform.

All four channels are synchronized and controlled together with Time and Power settings. The time range is 0.5 to 9.5 hours and is increased or decreased in half (0.5) hour increments. When the time displays 0.0, warming stops for all four (4) channels.

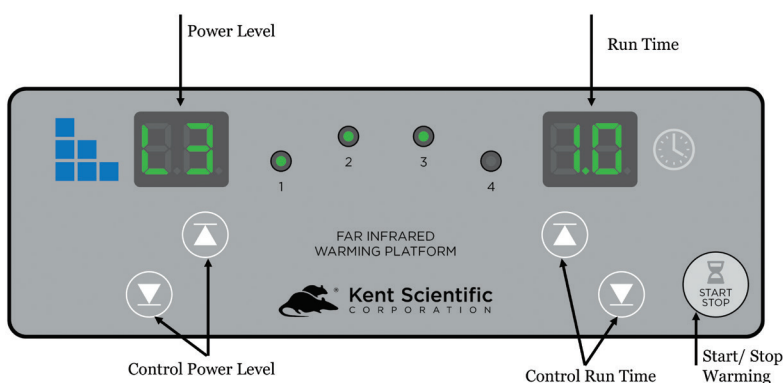

Temperature

1. Plug the power supply into the standard wall outlet. The unit will automatically power on.
2. Press the Up and Down buttons to select a warming level.

Tip: Set your platform to Level 3 then adjust as needed. There are many factors that affect the platform operating temperature. Use the included far-infrared thermometer and monitor animals' tail temperatures during use to ensure an acceptable temperature range.

3. Pre-warm the CODA[®] Animal Holders by placing the empty holders on the warming platform and cover with the CODA[®] Warming Cover.
4. Start the warming platform by pressing the START/STOP button on the platform.
5. Allow empty holders to pre-warm while configuring software settings.